


LIDT venting procedure:

- Turn off the flow bench

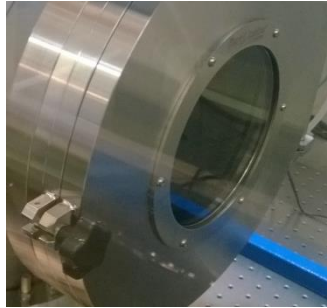


- Relies the lock on the chamber door

- Close the valve at the roughing pump



- Turn off the roughing pump (black switch on the back)
- Turn off the turbo pump by pressing star/stop button on the gray controller
- Wait until the turbo stops about half speed (you will hear loud noise)

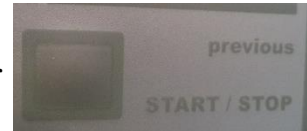

- Open the N tank with pressure set up to no more than 5 lb/inch²



- Open the valve on the gas line



- **VERY SLOWLY** open the up-to-air valve at the chamber



- Vent the chamber up the atmosphere by monitoring the pressure displayed by the gauges
- Immediately close the N tank, the valve at the gas line and the up-to-air valve at the chamber.

